



# Trigger Operated Eye, Face and Body Shower - Wall Mounting

Stock No. 763-486

A trigger operated eye, face and body shower incorporating a braided hose (1.5m long) and a Ballofix ball valve. To install the unit simply mount the wall bracket at a height of approximately 1-4 metres above floor level using the fixing screw and wall plug provided. A suitable drill for the purpose would be a 6mm masonry drill. Connect one end of the braided hose to the shower unit and the other end to the Ballofix ball valve. The water supply should then be connected and turned on.

Water inlet size 1/2" B.S.P. female. Minimum operating pressure 1.5 Bar G.

## Operating instructions

This trigger operated unit is essentially for use in flushing minor chemical splashes from the eyes, face, hands and other parts of the body.

The spray nozzle is fitted with a rubber cup to protect the user from further accidental injury. The nozzle is also protected against the ingress of dust and dirt by a dust cover, which is automatically removed by the initial jet of water. To operate the trigger operated unit simply remove the unit from the wall bracket, aim the nozzle at the contaminated area and actuate the hand trigger. For chemical in the eye, flush the open eye at once; continue for at least 5 to 10 minutes and, in any case of doubt, even longer. **If the contaminant is more than minimal seek medical advice as soon as possible.** The height of the spray can be regulated to suit the water supply pressure by adjusting the Ballofix ball valve using the Allen key provided.

## Maintenance

The unit should be cleaned regularly and inspected for any obvious damage. To clean first remove the protective rubber cap and dust cover from the main body, exposing the spray nozzle which should then also be unscrewed. The cap, dust cover and spray nozzle should be cleaned thoroughly. No sealants or solvents should be used. To refit, reverse the above operation.

Figure 1.

